




Automated Power Saturation Software

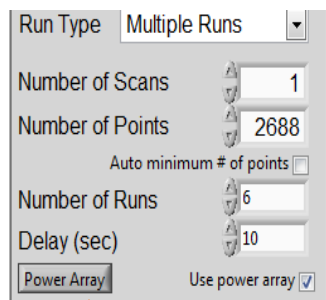

Power saturation data is key to obtaining good results for many ESR experiments. Active Spectrum's Micro ESR provides a simple automated way to acquire power data for any sample.

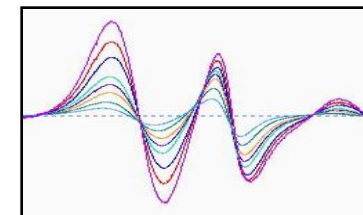


Signals from as asphaltene using an array of powers (above), Graph of amplitudes vs power (below).

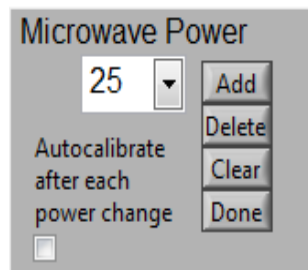


1. Just click the Power Array button when Multiple Runs has been selected.

2. The pop-up allows for simple setup of power array- Just select power from pull-down menu, and click Add.

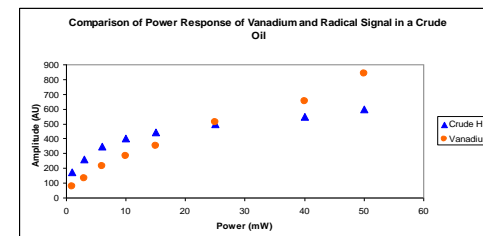

3. Choose as many powers as you like. Click Done when finished.

4. Option allows for tuning between runs.

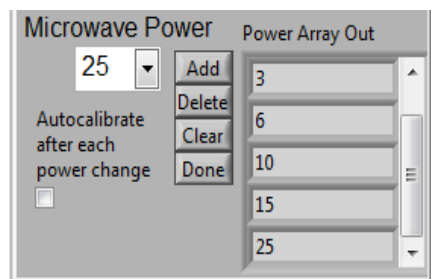

Power saturation data for various crude oils (below)

